

***Analyst™ Agilent Bug Fix Service Pack Release Notes***

*Version 1.3*

---

**General**

The Applied Biosystems/MDS SCIEX *Analyst™* Agilent Bug Fix service pack fixes the following bugs in Analyst 1.1:

- Difficulty in establishing communications with the Agilent devices from Analyst on Hardware Profile activation and communication problems with Agilent devices with newer version firmware (4.0x) and new board electronics.
- Entering -274 for the set point temperature in the Agilent 1100 column oven causes it to equilibrate forever.
- Does not turn off the temperature of the Agilent 1100 column oven when it starts up or when it goes into standby.

This service pack is intended for users who only want to fix the problems with Agilent devices in Analyst 1.1 without adding any new functionality. If you need control of the new Agilent devices (e.g., well plate autosampler, capillary LC pump, etc.) then please use the Agilent Upgrade service pack instead of this one.

The service pack requires 3 MB of disk space to install. This version of the service pack supercedes all previous versions and can be installed on top of any previous version. **Do not install this service pack on top of the Agilent Upgrade service pack.**

**Known Problems and Limitations**

Please report any problems, limitations and feature requests to Support via the Applied Biosystems/MDS SCIEX web site.

<http://www.appliedbiosystems.com/>

or via e-mail at:

[support@sciex.com](mailto:support@sciex.com)

The following are the known problems and limitations with this service pack:

- None

## **Installation Instructions**

**Before installing the service pack it is important to note that you must have administrator privileges for the computer you wish to install on. Contact your IT group if you are unsure what access rights you have.**

The installation of the Agilent Bug Fix is accomplished through the Agilent Bug Fix Installer. This program automatically places the required files in required folders. The Agilent Bug Fix cannot be uninstalled separately. It can only be uninstalled by uninstalling the *Analyst*<sup>TM</sup> Software. . The Agilent Bug Fix can be obtained from the Applied Biosystems website at: <http://www.appliedbiosystems.com/>

### **Step 1. Log on to the local system with administrator privileges.**

You must have administrator privileges on your local workstation in order to perform the installation.

### **Step 2. Installing the service pack software.**

Run the Agilent Bug Fix Installer, AgilentFix.exe.